


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: TRIGLYCERIDES (Olive Oil)

HPLC
Gradient

Fatty Acid Residues

OOL = Oleate/Oleate/Linolate
POL = Palmitate/Oleate/Linolate
OOO = Oleate/Oleate/Oleate
POO = Palmitate/Oleate/Oleate
SOO = Stearate/Oleate/Oleate

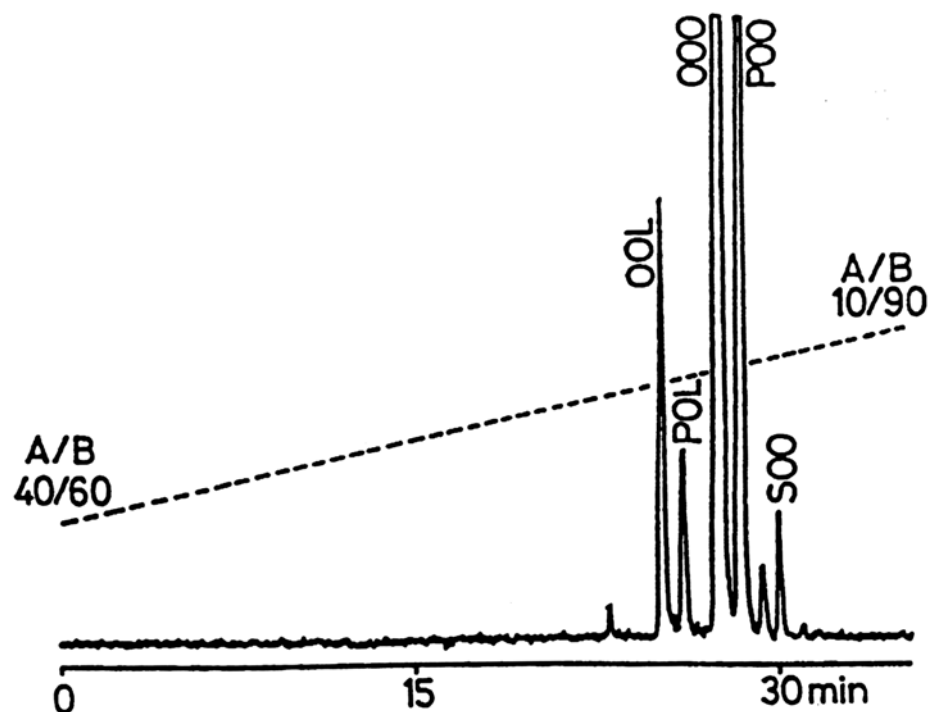

CONDITIONS

Column	Zorbax ODS - 7 μ m 250 x 9.4 mm
Injected Volume	1000 μ l
Eluent	A: Acetonitrile, B: CH ₂ H ₂
Flow Rate	1.5 ml/min
Temperature	30° C

Analyzed By : LCBA 45100 Orleans FRANCE [J. Chromatogr. 404 (1987) 95]