


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: ALKYL GLYCOSIDES

HPLC

- 1: Methyl α,β -D Glucopyranoside
- 2: Hexyl β -D Glucopyranoside
- 3: Octyl β -D Galactopyranoside
- 4: Octyl β -D Glucopyranoside
- 5: Octyl β -D Xylopyranoside
- 6: Decyl β -D Maltoside
- 7: Decyl β -D Glucopyranoside

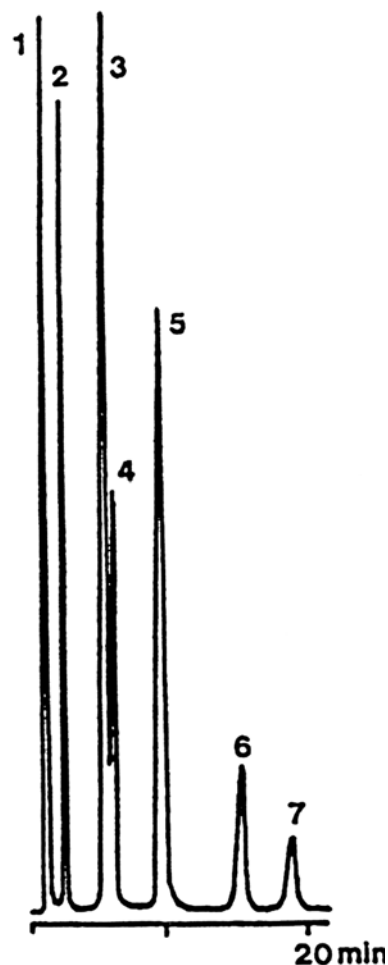

CONDITIONS

Column	Zorbax Phenyl - 7 μ m 15 x 4.6 mm
Injected Volume	20 μ l (Injected Quantity 8 μ g)
Eluent	Acetonitrile/Water (20/80)
Flow Rate	1.5 ml/min
Gas	Air
Pressure	2.5 Bars
Temperature	40° C

Analyzed By :LCBA 45100 Orleans FRANCE

S.E.D.E.R.E.

PARC VOLTA - BP 27

9, rue Parmentier

94141 Alfortville Cedex FRANCE

Application #26