


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: MALTODEXTRINS

HPLC
Gradient

Drug Excipient
Food Industry

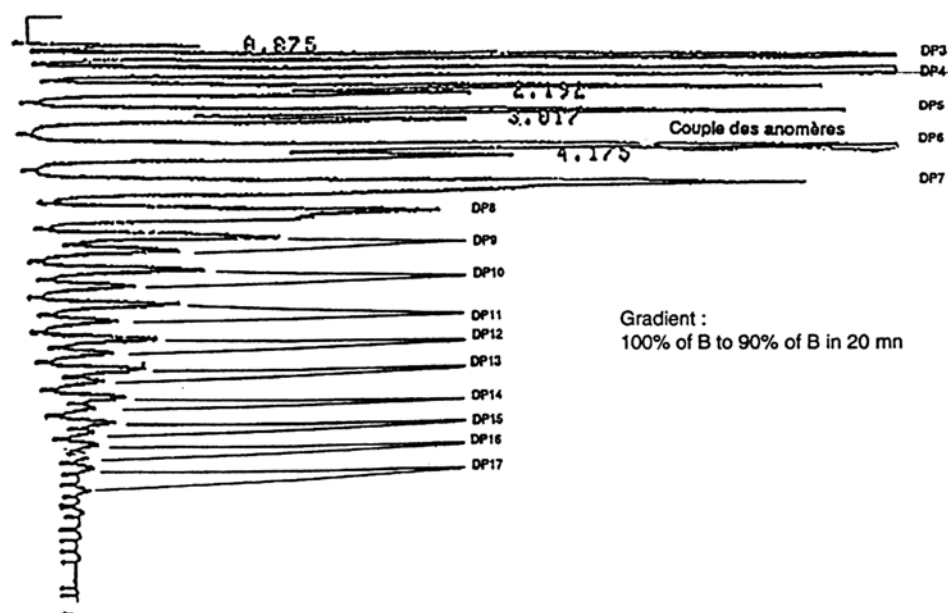

Gradient :
100% of B to 90% of B in 20 mn

Gradient: 100% to 90% of Water in 20 min

CONDITIONS

Column	Nucleosil C₁₈, 5 µm 15 x 4.6 mm
Injected Volume	20 µg, Injected Quantity: 1000 ppm
Eluent	A: Methanol, B: Water
Temperature	45° C