


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: MALTODEXTRINS

HPLC
Gradient

Drug Excipient
Food Industry

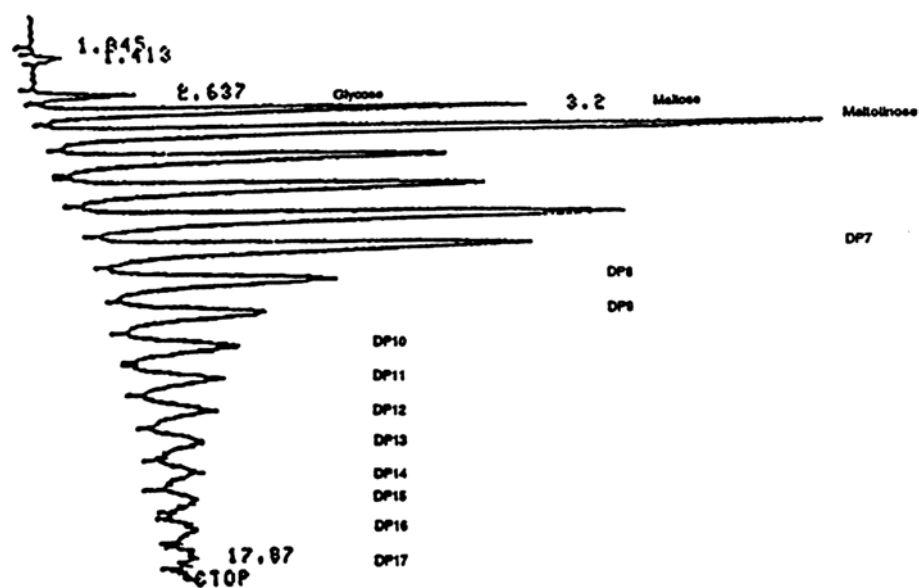

Gradient: 60 to 40 % Acetonitrile in 20 min

CONDITIONS

Column	Herbasil NH ₂ 250 x 4.6 mm
Injected Volume	20 µg, Injected Quantity 20 µg
Eluent	A: Acetonitrile, B: Water
Flow Rate	1.5 ml/min
Pressure	2.5 bars
Gas	Air
Temperature	40° C