


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: POLYETHYLENE GLYCOL 600

HPLC

Gradient

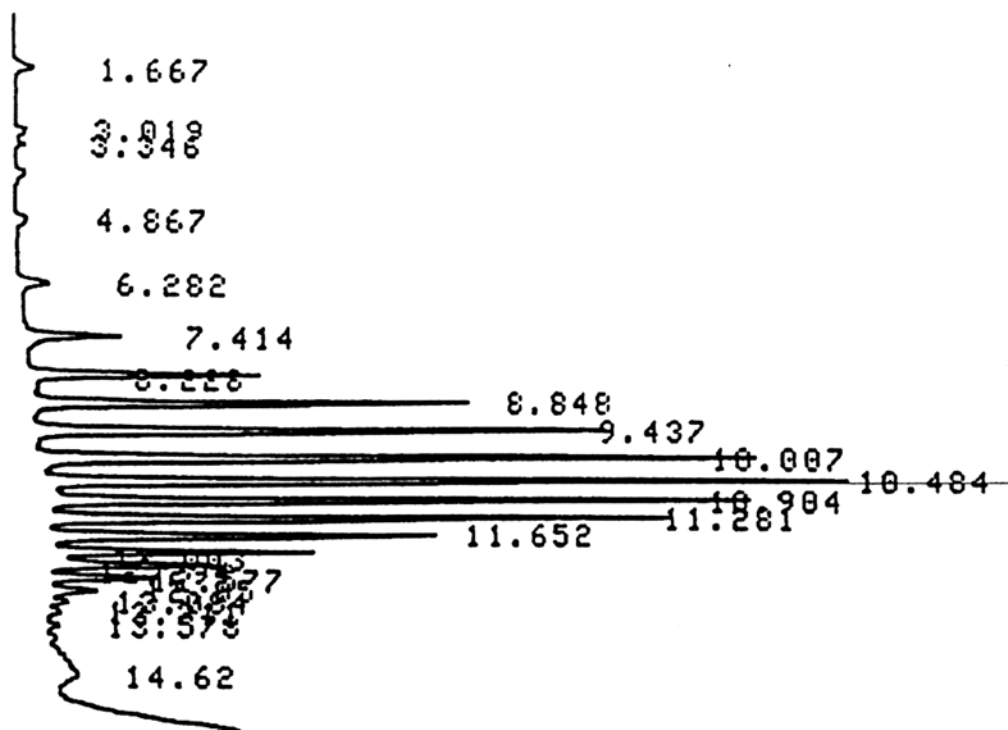

Gradient: 10% to 90% in 40 min

CONDITIONS

Column	<i>PRLPS 100 A, 5 μm</i> <i>250 x 4.6 mm</i>
Injected Volume	<i>20 μl, Injected Quantity 2 μg</i>
Eluent	<i>A: Acetonitrile</i> <i>B: Water</i>
Flow Rate	<i>1.3 ml/min</i>
Gas	<i>Air</i>
Pressure	<i>2.2 Bars</i>
Temperature	<i>40° C</i>

Analyzed by : LCBA 45100 Orleans FRANCE [J. Chromatogr. 505 (1990) 191]

S.E.D.E.R.E.

PARC VOLTA - BP 27

9, rue Parmentier

94141 Alfortville Cedex FRANCE

Application #52