


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: POLYETHYLENE GLYCOL 600

HPLC

Gradient

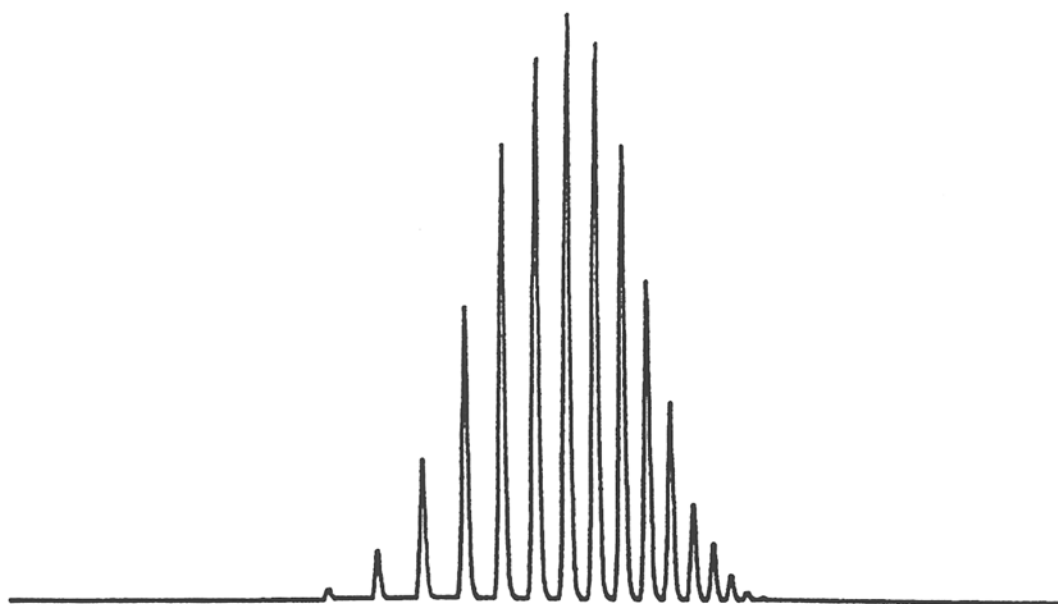

Gradient: 80% to 10% in 30 min
Hold at 10 % for 5 min
10 % to 80 % in 5 min

CONDITIONS

Column	<i>Nucleosil 120-7-C18</i> <i>250 x 4 mm</i>
Eluent	<i>A: Water/Methanol (95/5) [w/w/]</i> <i>B: Methanol/Water (95/5) [w/w/]</i>
Flow Rate	<i>1.0 ml/min</i>
Temperature	<i>36° C</i>