


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE: FATTY ACIDS (in Wine)

HPLC

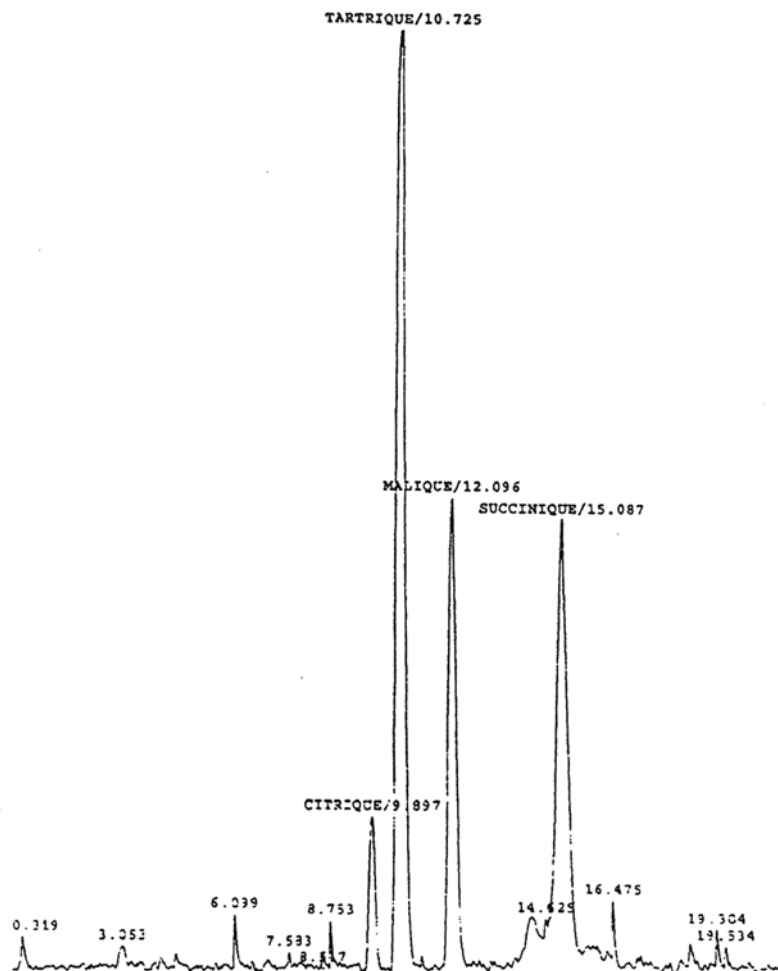

CONDITIONS

Column	CAR-H
Injected Volume	20 μ l
Eluent	Water + Trifluoroacetic Acid (pH = 2.3)
Flow Rate	1.0 ml/min
Pressure	2 bars
Gas	Air
Temperature	Column 40° C, Detector 40° C

Analyzed by: Touzart et Matignon, B.P. 52, 94403 Vitry Sur Seine FRANCE

S.E.D.E.R.E.
PARC VOLTA - BP 27
9, rue Parmentier
94141 Alfortville Cedex FRANCE

Application #71