


# SEDEX Low Temperature Evaporative Light Scattering Detector

**SAMPLE: ENANTIOMERS**

**HPLC**

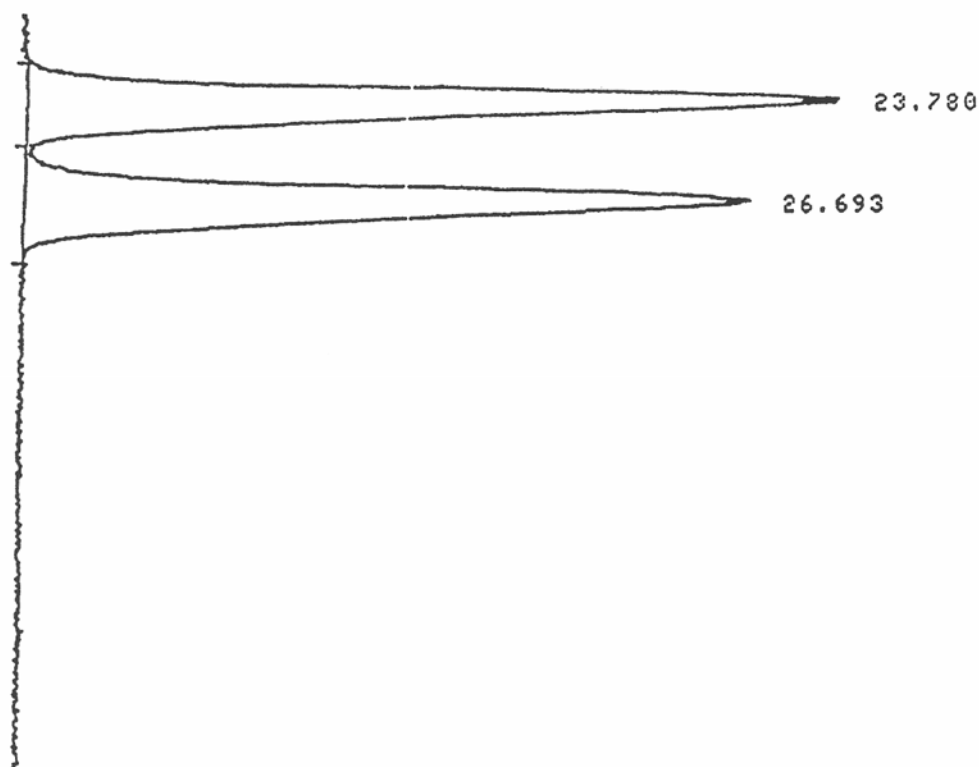

## CONDITIONS

|                 |                                           |
|-----------------|-------------------------------------------|
| Column          | <i>Chiracel OD</i><br><i>250 x 4.6 mm</i> |
| Injected Volume | <i>10 µl</i>                              |
| Eluent          | <i>Hexane/Ethanol (95/5)</i>              |
| Flow Rate       | <i>1.2 ml/min</i>                         |
| Pressure        | <i>2.0 bars</i>                           |
| Gas             | <i>Helium</i>                             |
| Temperature     | <i>40° C</i>                              |

Analyzed by: CE Saclay Molecules Marquees 91191 Gif Sur Yvette FRANCE

**S.E.D.E.R.E.**  
PARC VOLTA - BP 27  
9, rue Parmentier  
94141 Alfortville Cedex FRANCE

**Application #85**