


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE:

Micro HPLC

Gradient

- 1) Dimyristoylphosphatidyl glycerol, 1.2 µg/µl
- 2) Mannitol, 14 µg/µl
- 3) Dimyristoylphosphatidyl choline, 2.8 µg/µl

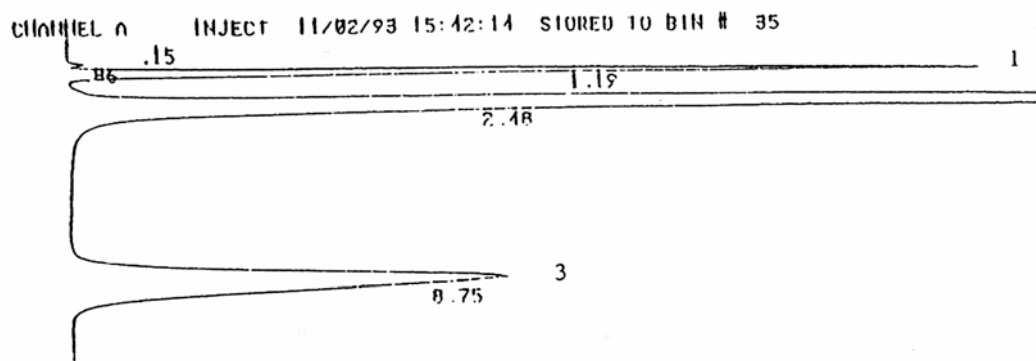

CONDITIONS

Column	Nucleosil Diol, 7 µm 250 x 4.0 mm
Injected sample	20 µl
Eluent	i-Propanol/Methanol/H ₂ O/NH ₄ OH (con) (57.5/36.5/5.9/0.1)
Flow Rate	1.2 ml/min
Pressure	2.2 bars
Gas	Nitrogen
Temperature	40° C

Analyzed By : Richard Scientific, Novato CA 94949 USA

S.E.D.E.R.E.

PARC VOLTA - BP 27

9, rue Parmentier

94141 Alfortville Cedex FRANCE

Application #95