


SEDEX Low Temperature Evaporative Light Scattering Detector

INORGANIC ANIONS

HPLC

Effect of drift tube temperature variation on sensitivity

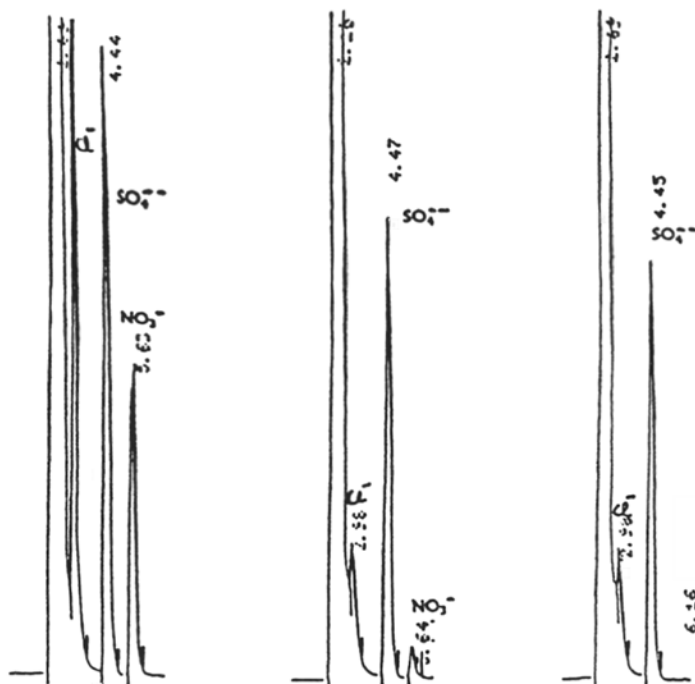

Stationnary phase :
HYPERCARB 7 μ m

Mobile phase :
HCOOH/NH₄OH

Sample :
PERRIER 20 μ l
Cl⁻ 22 ppm
NO₃⁻ 18 ppm
SO₄⁻ 33 ppm

CONDITIONS

Pressure	2 bars
Gas	Air
Gain	10
Temperature	45°C

Analyzed by : LCBA 45100 Orléans FRANCE

S.E.D.E.R.E.
PARC VOLTA - BP 27
9, rue Parmentier
94141 Alfortville Cedex FRANCE

Application # 108