


SEDEX Low Temperature Evaporative Light Scattering Detector

**SAMPLE: OLEINS,
TOCOPHEROLS CHOLESTEROL**

HPLC

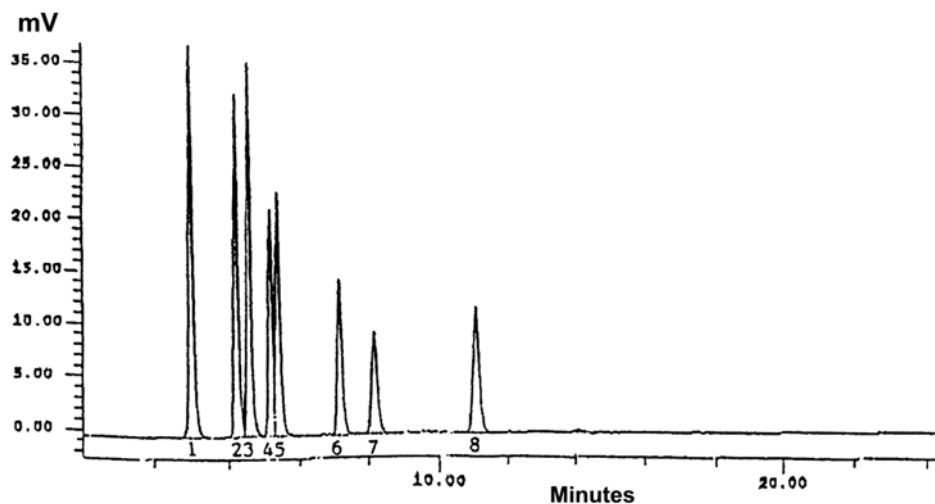

1. Triolein
2. α -Tocopherol
3. Oleic Acid
4. β -Tocopherol
5. γ -Tocopherol
6. δ -Tocopherol
7. Diolein
8. Cholesterol

CONDITIONS

Column	2 X Nova-Pak Si, 150 x 3.9 mm, 4 μ m
Injected volume	10 μ l 1,000 ppm ea.
mobile phase	Hexane/IPA (99/1)
Flow rate	1 ml/min
Pressure	2.2 bars
Gas	Nitrogen
Gain	6
Temperature	40°C

Analyzed by : Courtesy of Bunge Foods, Bradley, IL

S.E.D.E.R.E.
PARC VOLTA - BP 27
9, rue Parmentier
94141 Alfortville Cedex FRANCE

Application #116