


SEDEX Low Temperature Evaporative Light Scattering Detector

SAMPLE : PHENYLALANINE

HPLC

AMINO-ACID Direct Enantiomeric Separation

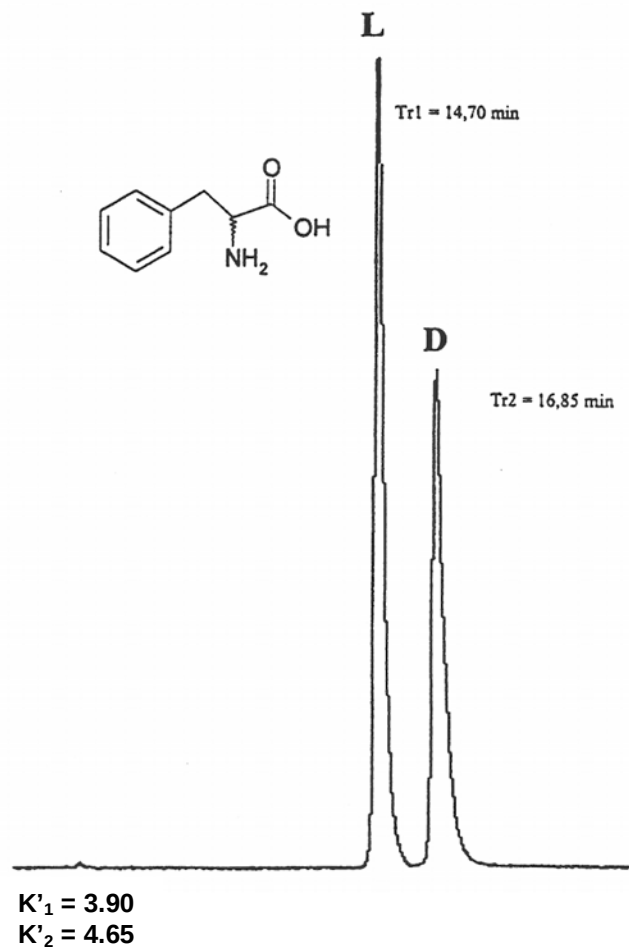

$\alpha = 1.19$
 $Rs = 2.80$

CONDITIONS

Column	Chirobiotic T 250 x 4.6 mm
Injected volume	5 μ l (1.5 ml)
mobile phase	Acetonitril / water 80/20
Flow rate	1 ml/min
Gas	Air
Gain	11
Temperature	25°C

Analyzed by : CEA SACLAY – Service des molécules marquées – 91191 GIF/YVETTE CEDEX FRANCE

S.E.D.E.R.E.

PARC VOLTA - BP 27

9, rue Parmentier

94141 Alfortville Cedex FRANCE

Application#127